

Glacios

45k Data Collection 0.872 Å/ubpix

LMM Settings

- Spot Size 6, 150 μm C2, LM 62 \times , C2 40.53%, Default Montage Settings

MMM Settings

- Spot Size 6, 50 μm C2, LM 380 \times , C2 70%

Low Dose Settings

- Spot size 6, 50 μm C2
- View: NanoProbe, SA 6700 \times , 79% C2
- Record/T/F: NanoProbe, SA 45k \times , 38.786% C2
- EPS 14, 0.877 Å/ubpix

Gain Ref Path

- C: drive>Program Data>Gatan>Reference Images
- File Format .dm4, most recent large file