

# Saving Ceta Camera Images in SerialEM

1. Start the serial EM server and Start SerialEM
2. Under Camera setup (Dark Green Control Panel) in serial EM, click Setup
3. Select Ceta camera, record image, gain Normalized, set binning and exposure time (This only needs to be done once)
4. Click OK at bottom of window
5. Take Record image, This is now your active image in buffer A
6. Click Save A or file Save active, specify file format, and file location within dosefractions
7. Once finished, the file will be open in the top left of SerialEM, hit file close and it will disappear
8. You are now ready to record and save your next image, all images will be transferred from your dose fractions folder to the corresponding folder on BioHPC.