


## FEG\_Registers

### For Krios ONLY:

LowDose -> FEG\_Registers -> select "300kV\_Dan\_..." -> click "Set"

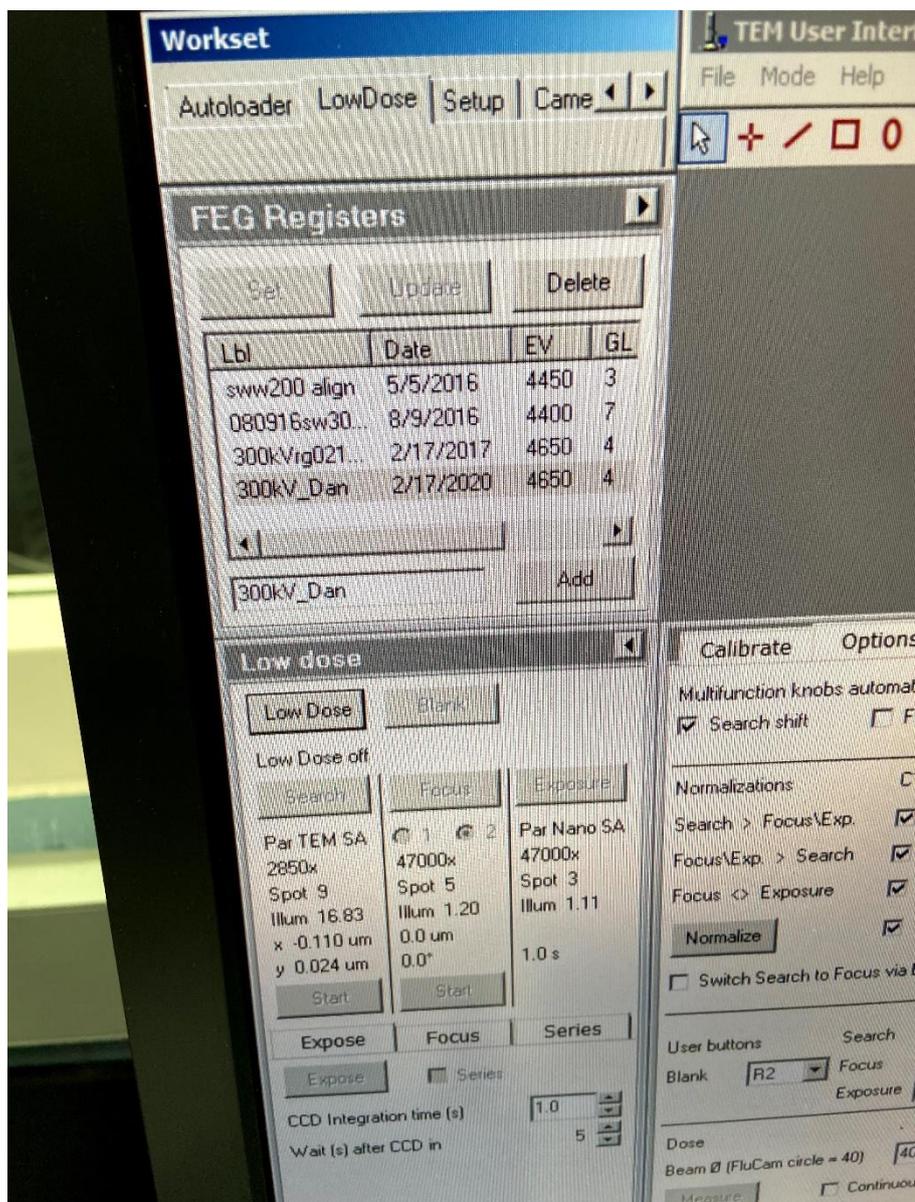

**Note:** the file navigator window on the right shows the location of the FEG\_Registers configuration files. To open it, click flap-out -> "File" -> "Open".

Users can save settings here as well.

## FEG\_Registers

### For Krios ONLY:

“Search” -> click “Alignments” -> flap-out -> click “File” -> select “300kV\_Dan\_...”

- Highlight all items on the bottom right corner window (green box) by click on top item, then hold the “SHIFT” key and click on the item at the bottom;
- click on the “left arrow” icon (blue arrow) so that same set of items show up on the window to the left;
- Click “Apply”.

